

Tips for Successful Cryopreservation

Technical Reference Guide

The practice of keeping a stock of frozen cultures should be a part of any cell culture laboratory. Cell lines or clones should be stored frozen to protect against genetic changes or loss of cultures due to contamination or incubator failure. Storage of cells in liquid nitrogen allows for prolonged storage, of up to 10 – 12 years. Cells can be stored in either the liquid nitrogen phase (-180°C to -195°C) or in the gas phase (-140°C to 145°C). The gas phase may be the better choice as the liquid can seep into the cryovial and cause the vial to “explode” upon thawing.

Cells to be frozen should be free of contamination and in the log phase of growth. The freezing medium should contain 5 – 10% cryopreservative and 5 – 20% serum. Commonly used cryopreservatives are glycerol or dimethyl sulfoxide (DMSO). Glycerol and DMSO lowers the freezing point. A slow cooling rate (1 – 2 degrees/minute) allows water to move out of the cell before it freezes. This protects against mechanical injury from ice crystal formation.

General Procedure

Adherent Cells

1. Remove cell culture medium.
2. Wash monolayer with calcium or magnesium-free PBS.
3. Remove PBS.
4. Add trypsin solution (just enough to cover monolayer).
5. Monitor the detachment process carefully to avoid cell damage.
6. When cells begin to round up, add complete medium containing serum or a trypsin-inhibitor solution.
7. Perform a cell count.
8. Cells will be frozen at 5×10^6 to 1×10^8 cells/ml (cryovials generally are 1 – 2 ml in size).
9. Based upon the number of vials to be prepared and cell number needed, centrifuge the appropriate volume of the cell suspension.
10. Aspirate off the growth medium.
11. Resuspend in freezing medium at 5×10^6 to 1×10^8 cells/ml.
12. Place vials in automatic freezing unit. If no such device is available, place vials on ice for about 1 – 2 hours, move to -20°C freezer for several hours, next move -70°C freezer for several hours, and finally into liquid nitrogen storage.

Suspension Cells

1. Perform a cell count.
2. Cells will be frozen at 5×10^6 to 1×10^8 cells/ml (cryovials generally are 1 – 2 ml in size).
3. Based upon the number of vials to be prepared and cell number needed, centrifuge the appropriate volume of the cell suspension.
4. Aspirate off the growth medium.
5. Resuspend in freezing medium at 5×10^6 to 1×10^8 cells/ml.
6. Place vials in automatic freezing unit. If no such device is available, place vials on ice for about 1 – 2 hours, move to -20°C freezer for several hours, next move -70°C freezer for several hours and finally into liquid nitrogen storage.

Thawing

1. Retrieve the cells from the freezer and quickly place it into a 37°C water bath and shake the vial. The cells must thaw quickly as the temperature from -50°C to 0°C has the greatest potential for cellular damage, so you want to have the temperature move quickly through that temperature range.
2. Remove cells from vial and slowly add growth medium until cells are at appropriate density for plating.
3. Plate in appropriate growth vessel.

In some cases, DMSO may be toxic to cells. For these, centrifuge cells for 2 minutes at 100 xg. Discard the freezing medium and resuspend the cells in fresh growth medium.

Contact Information

North America

Customer Service: 800 638 8174 (toll free)

order.us@lonza.com

Scientific Support: 800 521 0390 (toll free)

scientific.support@lonza.com

Europe

Customer Service: +32 87 321 611

order.europe@lonza.com

Scientific Support: +32 87 321 611

scientific.support.eu@lonza.com

International

Contact your local Lonza distributor

Customer Service: +1 301 898 7025

scientific.support@lonza.com

International Offices

Australia	+61 3 9550 0883
Belgium	+32 87 321 611
Brazil	+55 11 2069 8800
France	0800 91 19 81 (toll free)
Germany	0800 182 52 87 (toll free)
India	+91 40 4123 4000
Ireland	1 800 654 253 (toll free)
Italy	800 789 888 (toll free)
Japan	+81 3 6264 0660
Luxemburg	+32 87 321 611
Poland	+48 781 120 300
Singapore	+65 6521 4379
Spain	900 963 298 (toll free)
The Netherlands	0800 022 4525 (toll free)
United Kingdom	0808 234 97 88 (toll free)

Lonza Cologne GmbH – 50829 Cologne Germany

For research use only. Not for use in diagnostic procedures.
The Nucleofector™ Technology is covered by patent and/or patent pending rights owned by the Lonza Group Ltd or its affiliates.

Unless otherwise noted, all trademarks herein are marks of the Lonza Group or its affiliates. The information contained herein is believed to be correct and corresponds to the latest state of scientific and technical knowledge. However, no warranty is made, either expressed or implied, regarding its accuracy or the results to be obtained from the use of such information and no warranty is expressed or implied concerning the use of these products. The buyer assumes all risks of use and/or handling. Any user must make his own determination and satisfy himself that the products supplied by Lonza Group Ltd or its affiliates and the information and recommendations given by Lonza Group Ltd or its affiliates are (i) suitable for intended process or purpose, (ii) in compliance with environmental, health and safety regulations, and (iii) will not infringe any third party's intellectual property rights.