

Lonza Rockland, Inc. www.lonza.com biotechserv@lonza.com Tech Service: 800-521-0390 Customer Service: 800-638-8174 Document # 18924-1007-03 Rockland, ME 04841 USA

Formaldehyde Sample Buffer

For best results, use within 1 year from date of manufacture.

Recommended procedure:

- 1. Thaw RNA samples on ice.
- 2. Thaw sample buffer at room temp.
- 3. Vortex sample buffer to mix thoroughly.
- 4. Dilute RNA sample 1:1 with Formaldehyde Sample

Buffer.

- 5. Mix thoroughly.
- 6. Heat to 65°C for 5-15 minutes.
- 7. Cool for 1 minute on ice.
- 8. Load samples on gel.

Store: -20°C

For Research Use Only

© 2007 Lonza Rockland, Inc. All Rights reserved.