

Lonza Rockland, Inc. www.lonza.com biotechserv@lonza.com Tech Service: 800-521-0390 Customer Service: 800-638-8174 Document # 18148-0807-03 Rockland, ME 04841 USA

## Reliant® Gel System Reusable UV Transparent Tray

NOTE: remove tape from base of tray before use.

- 1. Remove Reliant® Precast Gel from disposable tray.
- 2. Place gel onto reusable tray.
- 3. Place this tray into the gel box oriented so that the open ends are toward the correct electrodes.
- 4. Electrophorese according to standard procedures.
- 5. Gel may be imaged in UV transparent tray.
- 6. Wash and store the tray for next use.

## **Ordering Information**

Catalog No.	Description	Size
54946	Portrait	Each
54945	Landscape	Each

For more information contact Technical Service at (800) 521-0390 or visit our website at www.lonza.com

All trademarks herein are marks of the Lonza Group or its affiliates.

© 2007 Lonza Rockland, Inc. All rights reserved.