



Lonza

# FlashGel<sup>®</sup> Cassette News

Simplified handling.

# Getting ready to load your samples has never been easier

## New seals on your FlashGel® Cassette.

There are still two seals on your FlashGel® Cassette: a larger white well seal with yellow tab and a smaller white vent hole seal with two black tips. This new seal concept allows the removal of well and vent seal in one step making cassette handling even more convenient.

### How does it work?

- Pull the yellow tab to remove well and vent seal
- Black tips are left behind on FlashGel® Cassette to partially cover vent openings

## Thank you.

For your trust in the FlashGel® System.

Lonza Rockland, Inc. – Rockland, ME 04841

Some components and technology of the FlashGel® System are sold under licensing agreements. The nucleic acid stain in this product is manufactured and sold under license from Molecular Probes, Inc. and the FlashGel® Cassette is sold under license from Invitrogen IP Holdings, Inc. and is for use only in research applications or quality control, and is covered by pending and issued patents.

For research use only. Not for use in diagnostic procedures. All trademarks belong to Lonza, registered in USA, EU or CH or to third party owners. The information contained herein is believed to be correct. No warranty is made, either expressed or implied. For more details: [www.lonza.com/legal](http://www.lonza.com/legal).