

# A Comparative Analysis of Absorbance-Based Endotoxin Testing Using Different Microplate Readers

Gerald Sando, David Santillan  
Lonza Walkersville, Inc., 8830 Biggs Ford Rd, Walkersville, MD 21793

## Purpose

- With Lonza’s planned discontinuation of sales of the ELx808™ Microplate Reader, commonly used for endotoxin testing, it is imperative for QC laboratories to find, evaluate, and implement a new reader for their normal operations. The aim of this study is to compare the performance of ELx808™ and Nebula® Absorbance Reader for their ability to measure absorbance-based endotoxin levels accurately and reliably
- Key Performance Indicators used for this study: accuracy, variability, and reaction time
- The results may be particularly useful for laboratories seeking to replace their current microplate readers or for those interested in optimizing their endotoxin testing protocols

## The Nebula® Absorbance Reader

- 96-well microplate reader
- Read capabilities:** Absorbance (monochromator)
- Light source:** Xenon Flash
- Detection:** Silicon photodiode
- Wavelength range:** 230 – 1000nm
- Optimized to work with WinKQCL® Endotoxin Detection and Analysis Software and traditional LAL assays such as the PYROGENT® 5000 Kinetic Turbidimetric Assay and the Kinetic-QCL® Kinetic Chromogenic Assay



## Results

Sample predictions


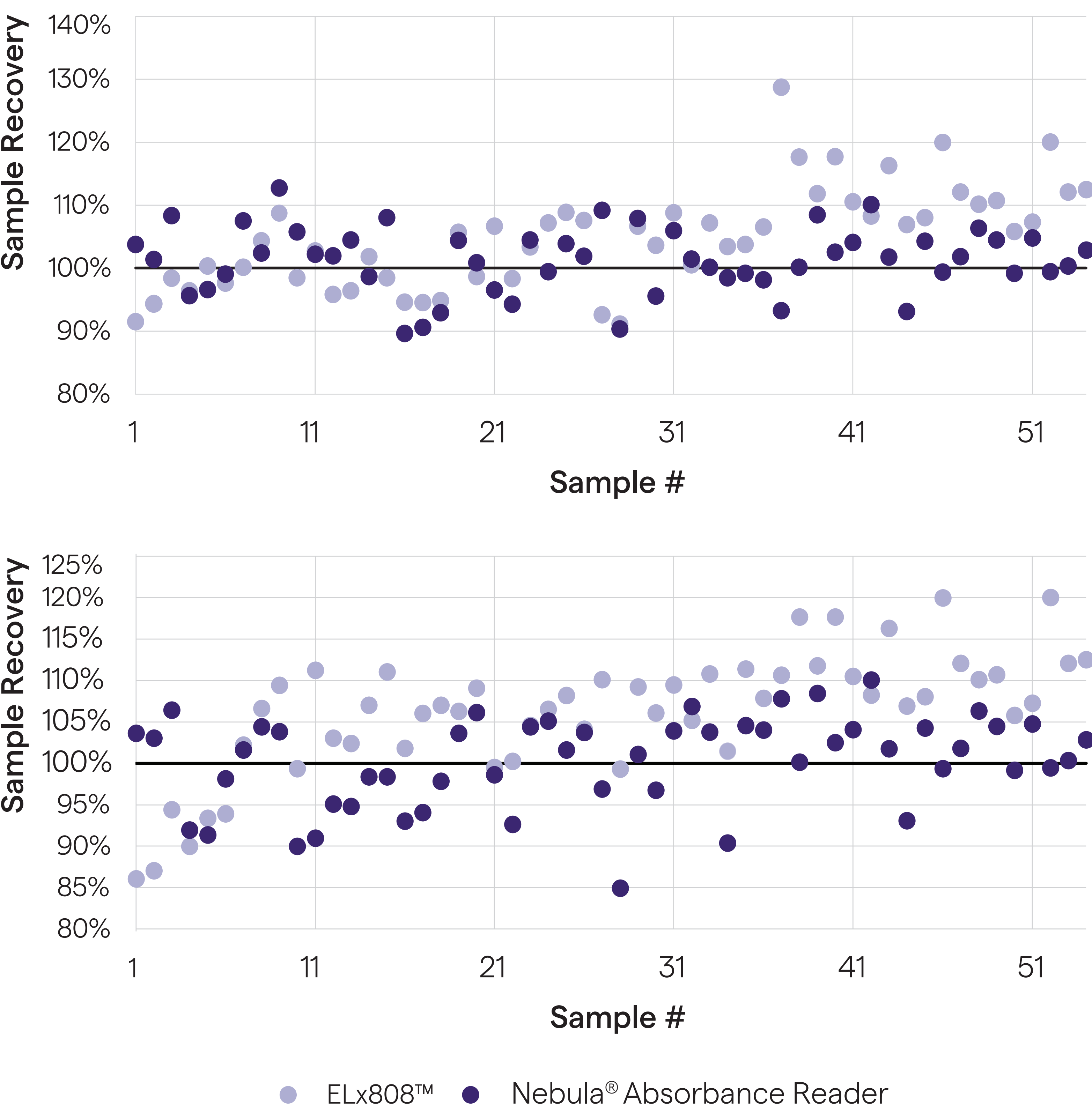


Figure 1. 54 samples for each chemistry, 18 at each concentration (0.025, 0.25, and 2.5 EU/mL for KQCL, 0.05, 0.5, and 5 EU/mL for Pyrogen® 5000)

	KQCL Sample Predictions		Pyrogen® 5000 Sample Predictions	
	Average	CV	Average	CV
ELx808™	105%	8.0%	106%	7.4%
Nebula® Absorbance Reader	101%	5.2%	100%	5.5%

## Uniformities

Kinetic-QCL® 0.5 EU/mL Uniformity Plate CVs

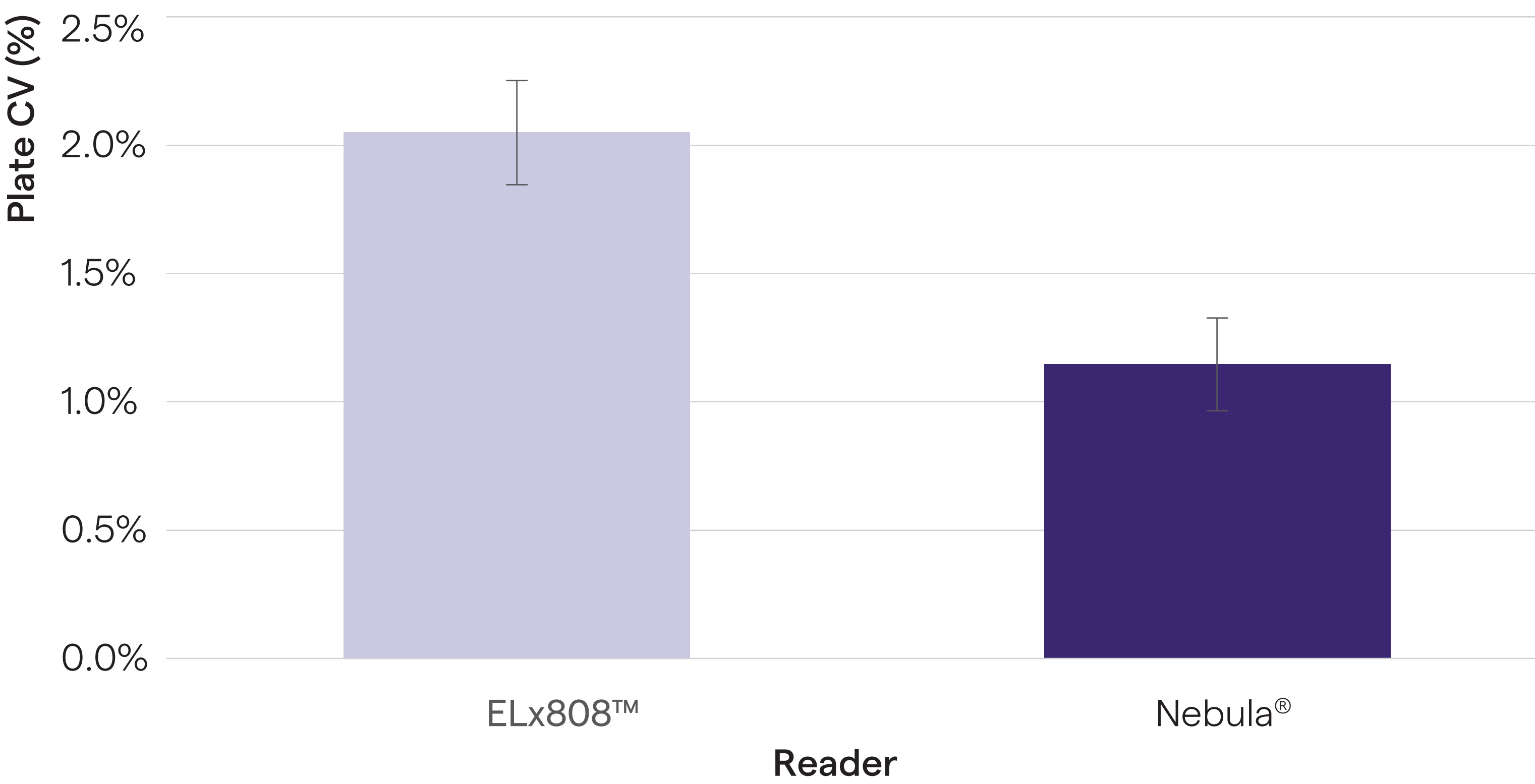

Pyrogen® 5000 1 EU/mL Unifirmity Plate CVs

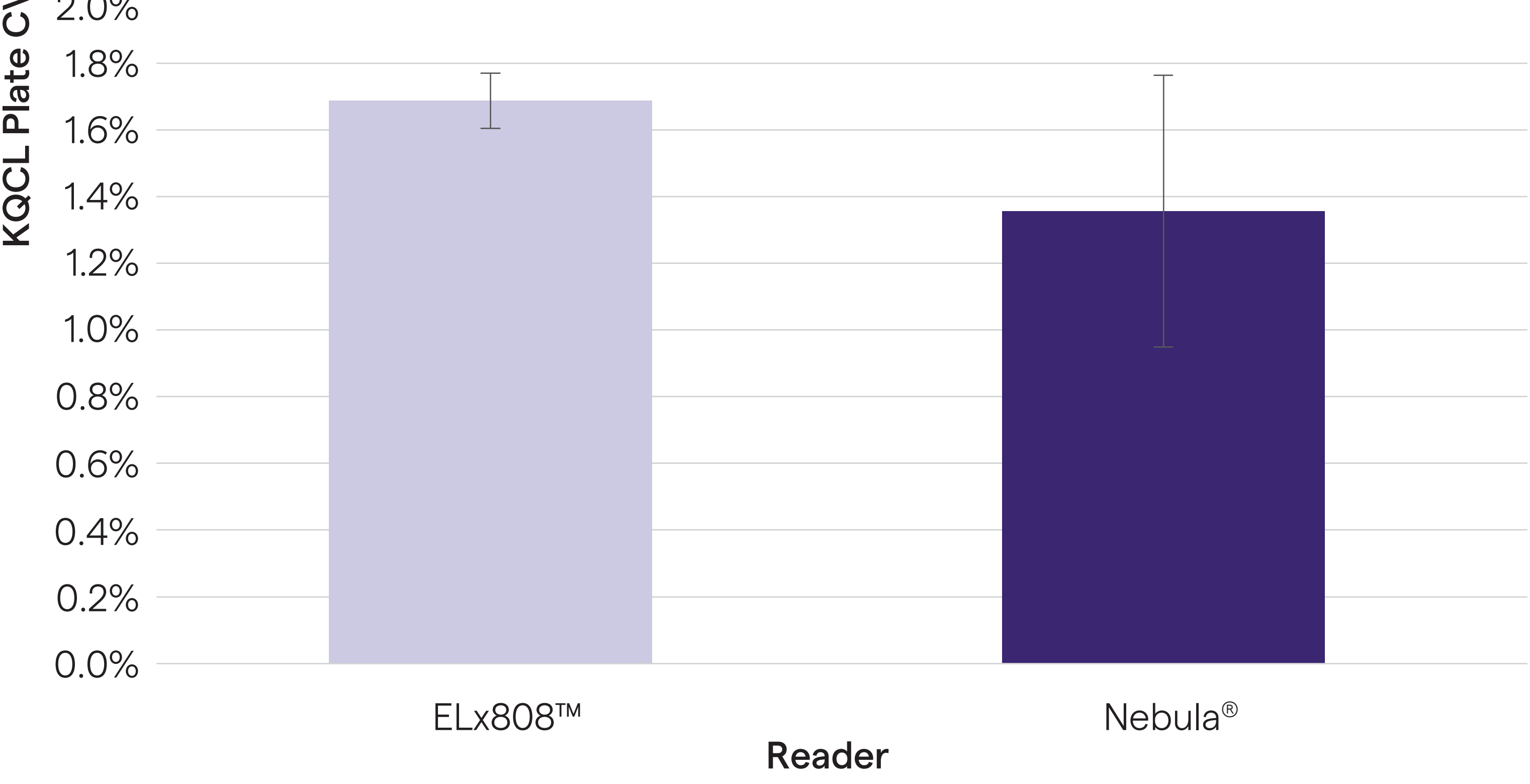



Figure 2. Uniformity testing

## Conclusion

- The data provided supports that the Nebula® Absorbance Reader is a suitable replacement for ELx808™ and it is fit for use for all of Lonza’s quantitative endotoxin detection absorbance assays
- Nebula® Absorbance Reader demonstrated comparable performance to ELx808™ with slight improvements in sample prediction accuracy and consistency

## Reaction Time

Kinetic-QCL® 0.005 EU/mL Reactions Times

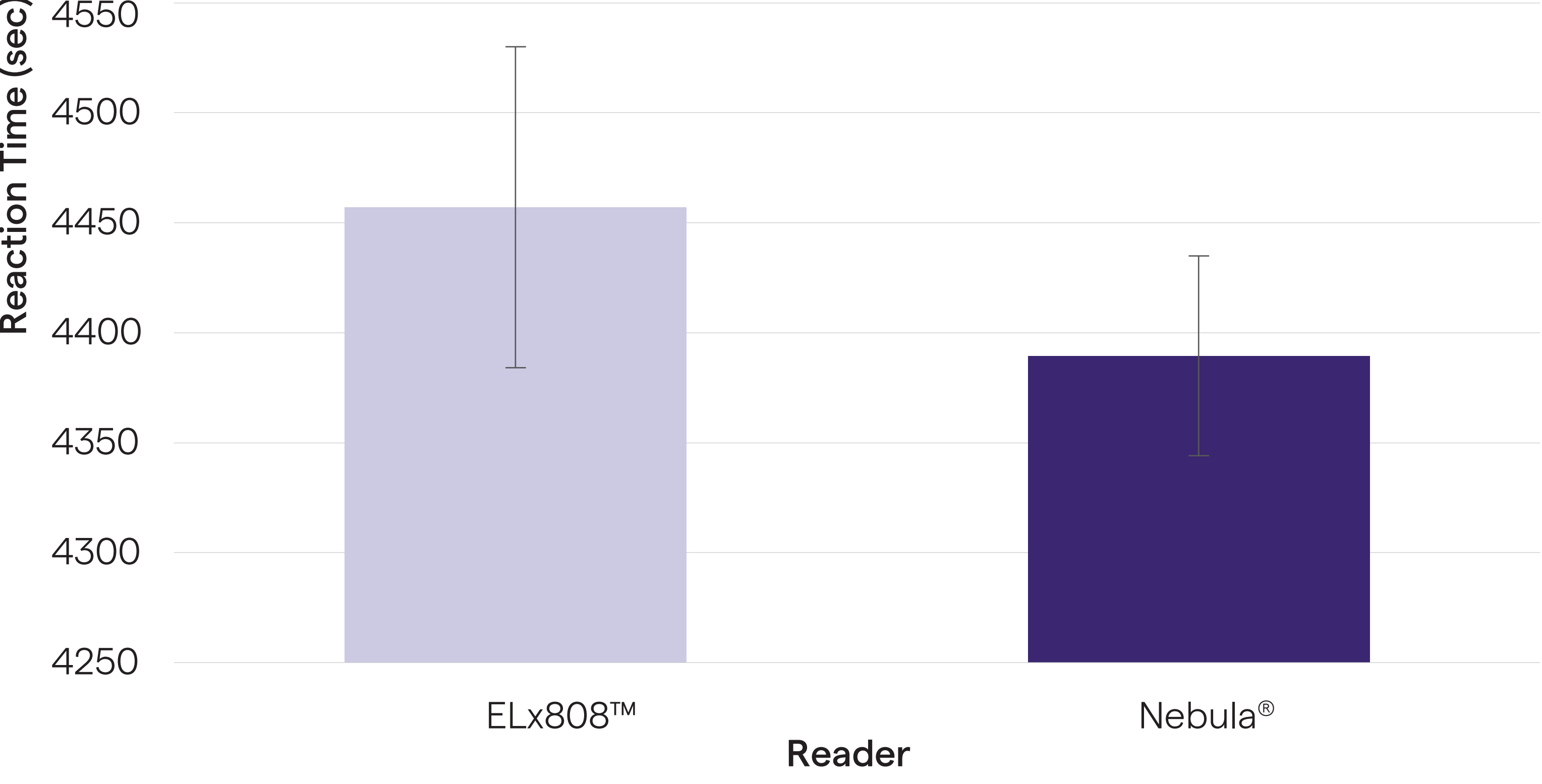

Pyrogen® 5000 0.01 EU/mL Reactions Times

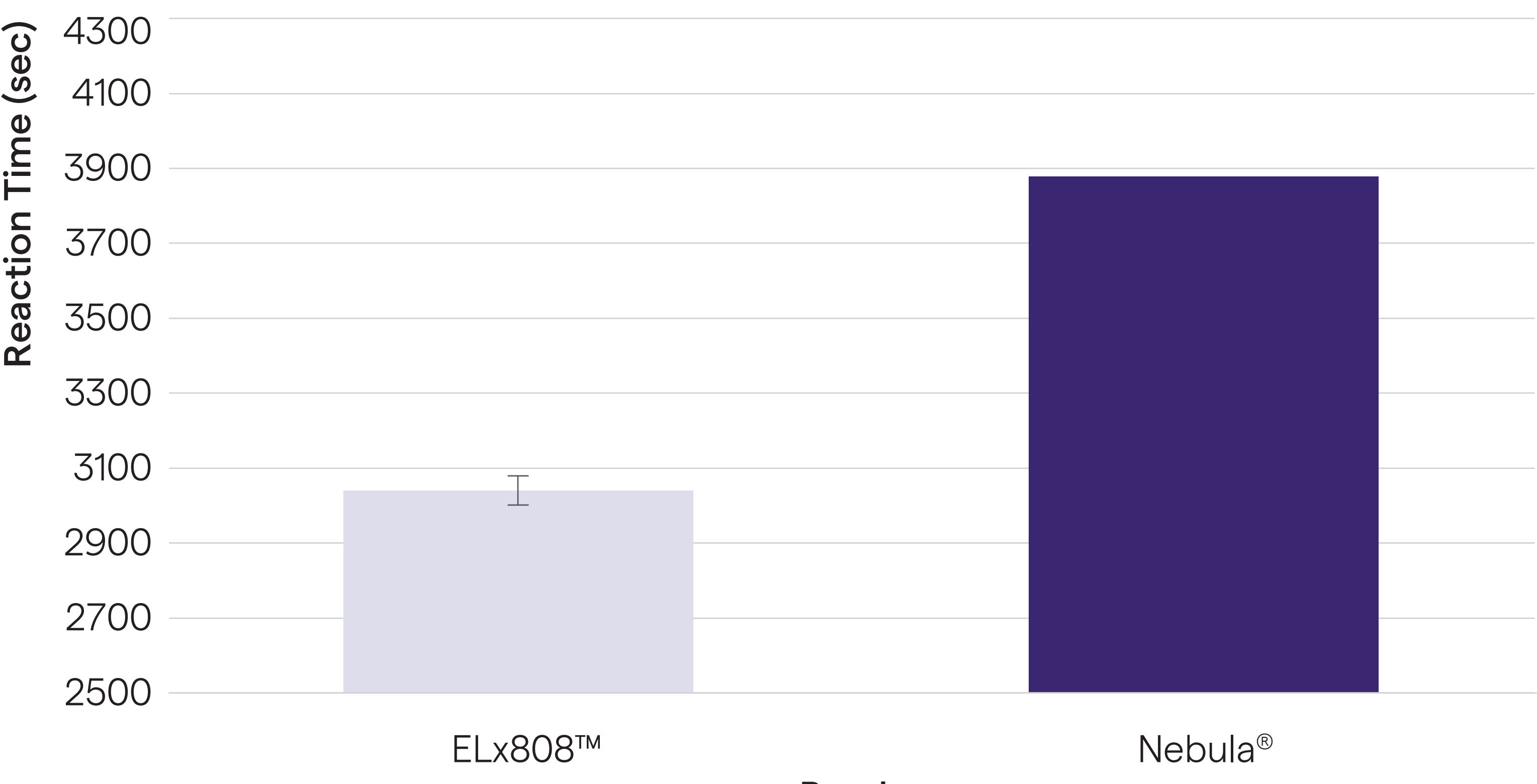


Figure 3. Reaction Time testing

- When using PYROGENT® 5000 Kinetic Turbidimetric Assay, the Nebula® Absorbance Reader demonstrated slower reaction times for the low standard (0.01 EU/mL)

All charts and figures were created using data from the Lonza’s 2023 Nebula Validation Study.  
All trademarks belong to Lonza, and are registered in USA, EU or CH or belong to third party owners and used only for informational purposes. The information contained herein is believed to be correct and corresponds to the latest state of scientific and technical knowledge. However, no warranty is made, either expressed or implied, regarding its accuracy or the results to be obtained from the use of such information and no warranty is expressed or implied concerning the use of these products. For more details: [www.lonza.com/legal](http://www.lonza.com/legal).  
©2023 Lonza, Inc. All rights reserved. RT-PO029 08/23