

4D-Nucleofector® LV Unit PRO

No commercial
use license fees



Efficient and scalable electroporation from research to GMP manufacturing

Experience the next generation 4D-Nucleofector® LV Unit PRO which comes with several new features and can be easily implemented into a GMP manufacturing process:

- Optimized cartridge design for improved performance for CRISPR-based knock-out/in or transposition
- Fixed volume cartridge with different filling volumes (0.5 mL, 1 mL, 1.5 mL and 2 mL) and new port for easy loading and clearing
- Flow-through cartridge with more robust sample processing
- Cartridge insertion support and visual acknowledgement

Transfection protocols can be established in smaller scale using the 4D-Nucleofector® X Unit and subsequently transferred to the LV Unit PRO with only a small re-optimization in 500 µL volume.

Benefit from

- Reliable, robust and efficient delivery of complex, clinically relevant cargos
- Easy scale up of the cell engineering process for up to 1 billion T cells
- Two vessel formats for closed transfection of medium or larger cell numbers
- Performance proven for resting and activated T cells
- Corresponding research-grade and GMP-grade consumables to transfer process development into GMP manufacturing
- 4D-Nucleofector® LogWare for operation under 21 CFR part11 compliant software
- Highly-skilled technical support
- Training support and IQOQ service

Find out more, visit:
lonza.com/transfection-scale-up

Robust scalability for different applications in activated T cells

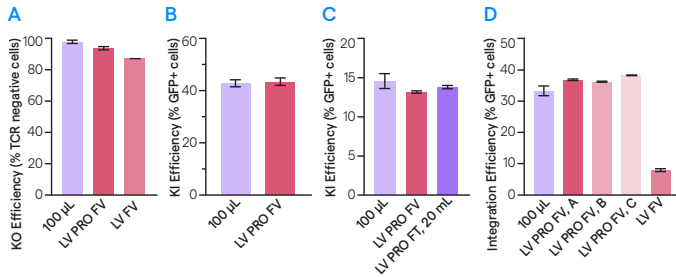



Figure 1. Comparable engineering efficiency for 100 µL Nucleocuvette® Vessel, fixed volume 2 mL Nucleocuvette® Cartridge PRO and flow-through LV Nucleocuvette® Cartridge PRO in activated T cells. (A) Knock-out of TRAC using Cas9 RNP (day 4, 2 donors, 4 technical replicates). (B) Knock-in of GFP into TRAC locus using Cas9 RNP and 3.5 kb dsDNA HDR template (day 4, 1 donor, 8 technical replicates). (C) Knock-in of GFP into RAB11A locus using Cas9 RNP and 3.5 kb dsDNA HDR template (day 4, 1 donor, 4 technical replicates). (D) Integration of GFP via piggyBac® transposase mRNA and transposon DNA with 3 different pulses (day 3, 1 donor, 4 technical replicates). The performance of the first-generation fixed volume 1 mL Nucleocuvette® Cartridge (LV FV) is included for comparison. Data are from representative donors. 100 µL = 100 µL Nucleocuvette® Vessel (100 µL), LV PRO FV = fixed volume 2 mL Nucleocuvette® Cartridge PRO; LV PRO FT = flow-through LV Nucleocuvette® Cartridge PRO. Days refer to post transfection.

Two vessel formats available in research or GMP quality



2 mL Nucleocuvette® Cartridge PRO

- Fixed volume samples of 0.5, 1, 1.5 or 2 mL
- Up to 2×10^8 cells
- Manual filling via optimized sterile injection port

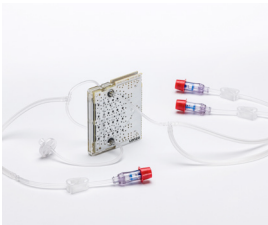

LV Nucleocuvette® Cartridge PRO

- Up to 20 mL processing volume
- Easily scalable to 1×10^9 cells
- Automatic filling via 4D-Nucleofector® LV Reservoirs or bags
- Weldable tubing

No commercial licensing fees

Cell therapy is already one of the most expensive treatments in history – why add additional costs? With Lonza's new 4D-Nucleofector® LV Unit PRO, you pay for the equipment and consumables. No commercial product use fees required.

A cost-efficient path to commercialization: help reduce your cost of goods and accelerate timelines without compromising quality.

Cat. no.	Description	Size
Instruments and accessories		
AAF-1003B	4D-Nucleofector® Core Unit	1 unit
AAF-1003X	4D-Nucleofector® X Unit	1 unit
AAF-1003L	4D-Nucleofector® LV Unit PRO Including 2 LV Reservoir Racks	1 unit
AAK-3001	4D-Nucleofector® LV Reservoir Rack	1 unit
SAAF-1002	4D-Nucleofector® LogWare (second generation)	1 unit
Research-grade consumables		
V4LP3-02250*	P3 Primary Cell Nucleofector® Solution LV Set	2.25 mL + 0.5 mL
V4LP3-22500*	P3 Primary Cell Nucleofector® Solution LV Set	22.5 mL + 5 mL
V4LN-8002	2 mL Nucleocuvette® Cartridge PRO	1 piece
V4LN-8020W	LV Nucleocuvette® Cartridge PRO, weldable tubing	1 piece
V4LR-1001W	4D-Nucleofector® LV Reservoirs, weldable tubing	2 pieces
GMP-grade consumables		
G4LP3-22500	TheraPEAK® P3 Nucleofector® Solution Set	22.5 mL + 5 mL
G4LP3-126000	TheraPEAK® P3 Nucleofector® Solution Set	126 mL + 28 mL
G4LN-8502	TheraPEAK® 2mL Nucleocuvette® Cartridge PRO - available end of 2025 -	5 pieces
G4LN-8520W	TheraPEAK® LV Nucleocuvette® Cartridge PRO, weldable tubing - available end of 2025 -	5 pieces
G4LR-1501W	TheraPEAK® 4D-Nucleofector® Reservoir, weldable tubing	5 pieces

* Corresponding SF Cell Line Solution Sets are available

Contact us

North America

Customer Service: +1 800 638 8174 (toll free)
 order.us@lonza.com
 Scientific Support: +1 800 521 0390 (toll free)
 scientific.support@lonza.com

Europe

Customer Service: +32 87 321 611
 order.europe@lonza.com
 Scientific Support: +49 221 99199 400
 scientific.support.eu@lonza.com

International

Contact your local Lonza distributor
 Customer Service: +1 301 898 7025
 scientific.support@lonza.com

Lonza Cologne GmbH – 50829 Cologne, Germany

Learn more.



Lonza Group Ltd. and its affiliates (collectively and individually, "Lonza") make efforts to include accurate and up-to-date information. However, Lonza makes no representations or warranties, express or implied, including as to accuracy or completeness of information. All trademarks belong to Lonza, and are registered in the USA, EU and/or CH, or used in common law, or belong to third-party owners and are used for only informational purposes. All third-party copyrights have been reproduced with permission from their owners. The user bears the sole responsibility for determining the existence of any third-party rights and obtaining any necessary licenses and approvals. For more information, including regarding legal disclaimers, Lonza's intellectual property rights, and how Lonza collects, uses and protects personal information: www.lonza.com/legal, <https://www.lonza.com/about-us/strategy/intellectual-property> and <http://www.lonza.com/privacy>.