




Top 10 tips for Mycoplasma free cell culture

Wear appropriate personal protective equipment (gloves, dedicated clean lab coat, safety glasses) and disinfect gloves

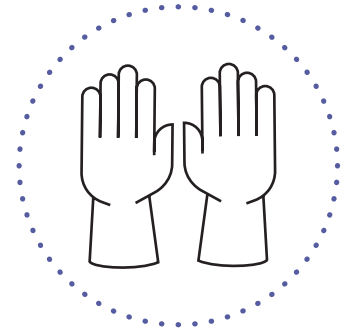

Wipe down work surfaces in safety cabinet before starting and in between cell lines

Avoid talking over your cells or sneezing

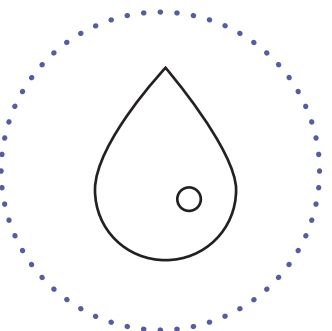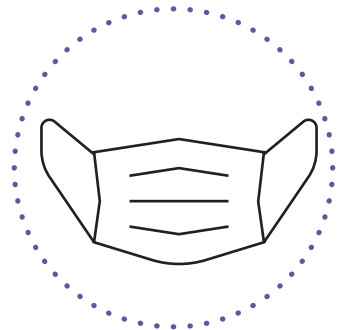



Avoid splashes, spills and aerosols

Use designated media bottles for each cell line




Clean incubator including water pan on a regular basis

Empty and clean water bath regularly

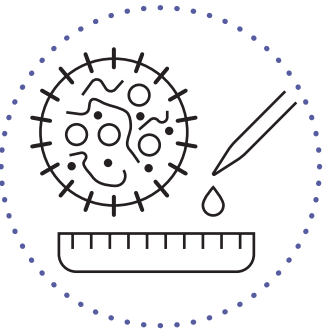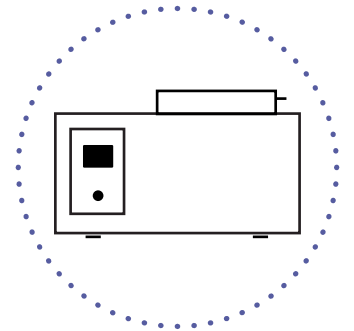

Test new cell cultures for mycoplasma contamination a few days after thawing

Quarantine all cell cultures coming into the lab until they have been tested for mycoplasma

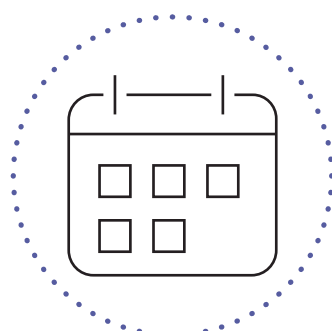

Don't forget: Perform regular mycoplasma testing every other week

More information :

www.lonza.com/mycoplasma

