

Lonza Walkersville, Inc. www.lonza.com scientific.support@lonza.com Scientific Support: 800-521-0390 Document # TS-CC-44-10 11/10 Walkersville, MD 21793-0127 USA © 2010 Lonza Walkersville, Inc.

Clonetics™ Bronchial Smooth Muscle Cell Systems

BSMC & DBMSC



Clonetics™ Bronchial Smooth Muscle Cell Systems contain Normal Human Bronchial Smooth Muscle Cells (BSMC) or Diseased Human Bronchial Smooth Muscle Cells (DBSMC) and optimized media for their growth. Each System can generate BSMC cultures for experimental applications in respiratory disorders including asthma, cystic fibrosis and respiratory distress syndrome. Clonetics™ Bronchial Smooth Muscle Cell Systems are convenient and easy to use, allowing the researcher to focus on results. There are three types of DBSMC available – cells isolated from donors with cystic fibrosis, asthma, and COPD. Cryopreserved BSMC and DBSMC are shipped in third passage. Proliferating BSMC and DBSMC are shipped in fourth or fifth passage.

Clonetics[™] Cells, Medium and Reagents are quality tested together and guaranteed to give optimum performance as a complete Cell System.

Cell System Components (Need to be purchased separately)

- One Bronchial Smooth Muscle Cell Product Normal or Diseased (Cryopreserved or Proliferating)
- One Smooth Muscle Cell Medium BulletKit™ -500 ml
 - Clonetics™ SmGM™-2 BulletKit™ (CC-3182) contains one 500 ml bottle of Smooth Muscle Cell Basal Medium and the following growth supplements: hEGF, 0.5 ml; Insulin, 0.5 ml; hFGF-B, 1ml; FBS, 25 ml; GA-1000, 0.5 ml.
- One ReagentPack™ (CC-5034) Containing:
 Trypsin/EDTA 100 ml
 Trypsin Neutralizing Solution 100 ml
 HEPES Buffered Saline Solution 100 ml

Characterization of Cells

Routine characterization of BSMC and DBMSC includes immunofluorescent staining. Cells stain positive for smooth muscle α -actin and negative for

von Willebrand (Factor VIII) antigen after differentiation.

Performance

Recommended seeding density for subculturing	3,500 cells/cm ²
Typical time from subculture to confluent monolayer	6 - 10 days
BSMC additional population doublings guaranteed using Clonetics™ System	15
DBSMC additional population doublings guaranteed using Clonetics™ System	Tested through 3 passages for information only

Quality Control

HIV-1, Hepatitis B and Hepatitis C are not detected for all donors and/or cell lots. All cells are performance assayed and test negative mycoplasma, bacteria, yeast and fungi. Cell viability, morphology and proliferative capacity are measured after recovery from cryopreservation. Clonetics™ Media are formulated for optimal growth of specific types of normal human cells. A Certificates of Analysis (COA) for each lot of cryopreserved cells is shipped with each order. COAs for all other products are available upon request.

Ordering Information

Cryopreserved Cells

CC-2576	BSMC	≥ 500,000 cells
00196980	DBSMC-CF (cystic fibrosis)	≥ 500,000 cells
00194850	DBSMC – AS (asthma)	≥ 500,000 cells
00195274	DBSMC-COPD	≥ 500,000 cells

Proliferating formats are available. Refer to the Lonza website or contact Scientific Support for details.

Lonza

CC-3182	SmGM [™] -2 BulletKit [™] , SmBM [™] plus SingleQuots [™] of Growth Supplements	500 ml
CC-3181	SmBM™, Smooth Muscle Cell Basal Medium	500 ml
CC-4149	SmGM [™] -2 SingleQuots [™] , Formulates SmBM [™] to SmGM [™] -2	
CC-5034	ReagentPack™	
	Trypsin Neutralizing Solution	100 ml
	Trypsin/EDTA Solution	100 ml
	HEPES Buffered Saline Solution	100 ml

When placing an order or for technical service, please refer to the product numbers and descriptions listed above. For a complete listing of all CloneticsTM Products, refer to the Lonza website or the current Lonza catalog. To obtain a catalog, additional information or technical service you may contact Lonza by web, e-mail, telephone, fax or mail.

Product Warranty

CULTURES HAVE A FINITE LIFESPAN *IN VITRO*. Lonza guarantees the performance of its cells only if Clonetics™ Media and Reagents are used exclusively, and the recommend protocols are followed. The performance of cells is not guaranteed if any modifications are made to the complete Cell System. Cryopreserved BSMC and DBSMC are assured to be viable and functional when thawed and maintained properly.

THESE PRODUCTS ARE FOR RESEARCH USE ONLY. Not approved for human or veterinary use, for application to humans or animals, or for use in clinical or *in vitro* procedures.

WARNING: CLONETICS™ AND POIETICS™ PRODUCTS CONTAIN HUMAN SOURCE MATERIAL, TREAT AS POTENTIALLY INFECTIOUS. Each donor is tested and found non-reactive by an FDA approved method for the presence of HIV-I, Hepatitis B Virus and Hepatitis C Virus. Where donor testing is not possible, cell products are tested for the presence of viral nucleic acid from HIV, Hepatitis B Virus, and Hepatitis C Virus. Testing can not offer complete assurance that HIV-1, Hepatitis B Virus, and Hepatitis C Virus are absent. All human sourced products should be handled at the Biological Safety Level 2 to minimize exposure of potentially infectious products, as recommended in the CDC-NIH Manual, Biosafety in Microbiological and Biomedical Laboratories, 5th Edition. If you require further information, please contact your site Safety Officer or Scientific Support.